Synchro set

Protocollo di utilizzo
Protocole d'utilisation
Benutzungsprotokoll
Protocol of use
Protocolo de utilización
Protocolo de utilização
Πρωτόκολλο χρήσης

- PER COLTURADI CELLULE
- POUR LA CULTURE DES CELLULES
- FÜR ZELLKULTUREN
- FOR CELL CULTURE
- PARA CULTIVO DE CÉLULAS
- PARA CULTURA DE CÉLULAS
- ΓΙΑ ΚΑΛΛΙΕΡΓΕΙΑ ΚΥΤΤΑΡΩΝ

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Synchronization kit for peripheral blood and marrow cells.

code EK AMTS 008 50 tests

Supplied material

- 4 vials x 1.5 ml of solution A
- 4 vials x 1.5 ml of solution B

Conservation and stability

Synchroset is shipped at +4°C. Once the product is received, it must be stored at -20 ± 2°C. The shelf life of the product is 12 months. Synchroset can be defrosted and frozen without altering its functionality.

English

Protocol of use

Culture synchronization

1. After 48-72 hours of culture, add 20 µl of solution A for ml of suspension fluid.
2. Incubate overnight.
3. Add 20 µl of solution B for ml of suspension fluid.
4. Incubate for 5 hours (it is not necessary to wash the cells).
5. Add 4 µl of Colcemid (10 µg/ml) for ml of suspension fluid.
6. Incubate at 37 ± 2°C for about 1 hour. Incubation may be reduced to 15 minutes to obtain a high number of prometaphases.

The chromosome mixes, after cell synchronization, require longer treatment times with trypsin to obtain a suitable GTG or RHG banding technique.

NB: for high quality results, it is strongly suggested to use Synchroset in combination with EuroClone media optimized for cell synchronization. Chromosome Synchro P and M are kits specifically realized in order to attain an efficient synchronization of the cell cycle, for culture and following analysis of samples of peripheral blood and marrow.

Cell synchronization which can be attained with Synchroset together with accurate formulation of properly studied media, guarantee an excellent mitotic index and high banding resolution (up to 550 bands for kit) thanks to a synergic effect of the two components.

Code	Description	Format	Use
EKAMTSY20	Chromosome Synchro P	20 culture tubes	Kit for culture and synchronization of peripheral blood cells
EKAMTSY50		50 culture tubes	
EKAMTBSY100.2		2 bottles of 100 ml	
EKAMTBSY100.5		5 bottles of 100 ml	
EKAMTBSY500		1 bottle of 500 ml	
EKAMTSY20M	Chromosome Synchro M	20 culture tubes	Kit for culture and synchronization of marrow cells
EKAMTSY50M		50 culture tubes	
EKAMTBSY100.2M		2 bottles of 100 ml	
EKAMTBSY100.5M		5 bottles of 100 ml	
EKAMTBSY500M		1 bottle of 500 ml	

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EuroClone S.p.A.